# Safeguarding Hygiene Where it Matters Most





# **Lumitester** ™ **Smart**

Hygiene monitoring in 10 seconds



**HYGIENE MONITORING IN** THREE EASY STEPS



**VISUALIZE THE DATA VIA APP** 



**ACCESS FROM ANYWHERE WITH CLOUD STORAGE** 

+ LuciPac TM With A3 Detection

ATP + ADP + AMP

More Reliable and Sensitive than ATP Tests A3 Finds What Others Miss



www.hyserve.com



info@hyserve.com



+49(0)89 3703 1223







# MAKE HYGIENE MONITORING A HABIT FOR EVERYONE

With the App, anyone can check cleanliness easily



Easy-to-use, results within 10 seconds



#### Swab

Swab the sample on a 10 x 10cm area, and insert the swab into main body



### Measure

Shake well and insert into the Lumitester





**Expiry** 

## Analyze

Measurements (RLU) displayed within 10 seconds



App to automatically analyse data

Recorded data can be displayed in time-series by registering with the Lumitester Smart App. Inspection pass rates are automatically graphed and data can be visualized. Employees will be more conscious of hygiene, and will maintain a high standard of environmental sanitation, which will help build greater trust in your business.







Visualize overall inspection scores through graphs



Cloud data management to combine data from multiple facilities

Data will be stored in the cloud and can be accessed from anywhere in the world. This will allow fast detection and response to any problems that may arise.



## **Lumitester** <sup>™</sup> **Smart** Product code: 1902653

Measurement time	10 seconds
Detection	ATP + ADP + AMP
Detection method	Luciferase-Enzyme Assay
Unit of Meassurement	Relative Light Units (RLU)
Included Accessories	2 alkaline batteries (AA), Cleaning brush, USB cable, Strap, Quick Manual

- Use for hygiene monitoring and other applications where detection of organic residues needs to be performed.
- Lucipac A3 and Lumitester Smart should not be used for counting general living bacteria or detecting specific pathogens.

LuciPac™ A3 Surface LuciPac™ A3 Water	Product code: 1702671 Product code: 1702672	
Storage Condition	2-8°C (do not freeze)	
25°C	14 days (unopened)	
30°C	5 days (unopened)	

15 months after manufacturing date



HyServe GmbH & Co KG. Ammerthalstraße 7 Tel. +49 (0) 89 3703 1223

info@hyserve.com

D- 85551 Kirchheim/München Fax +49 (0) 89 3703 1225 www.hyserve.com